



Chocolate Analyzer

EQUIPMENT SERIES



- Cocoa and Chocolate Ingredients
- Fat, Moisture, Sugar, Fatty Acids and Other Parameters in Under 30 Seconds
- Ready to go Pre-Calibrated with Custom Calibrations for Chocolate Industry



Process and Quality Control

Fast, accurate analysis is critical for both cocoa processors and chocolate manufacturers to optimize their process and insure quality. Optimizing the process will save the plant time and money and provide very rapid payback.

Near Infrared (NIR) analysis is a proven technique designed to provide fast, accurate, and reliable results for the cocoa industry.

BENEFITS:

- Rapid payback
- At-line ready
- Simple to use
- Flexible sample handling
- Easy data integration
- Low maintenance
- ISO 12099 Compliant
- Guaranteed calibrations

Monitoring incoming cocoa beans and other ingredients for quality

Analyze nibs, cocoa liquor and cocoa powder in process for quality and consistency

Verify fat, cocoa and sugar content of final chocolate products



SpectraStar™

The Unity Chocolate Analyzer is based on the SpectraStar XT, a state of the art hi-performance NIR analyzer optimized for the analysis of ground and unground natural products. The SpectraStar XT contains TAS, a revolutionary technology which monitors the instrument status and precisely calibrates the instrument to first principles. The SpectraStar XT delivers reliable operation and ultimate performance.

The Unity Scientific SpectraStar™ RTW (Rotating Top Window) system gives the ultimate in sampling flexibility and allows the customer to make rapid measurements of all important processing steps.

Plug and Go Calibrations

The Unity Chocolate Analyzer has ready to use calibrations for every step of the chocolate manufacturing process, from incoming cocoa beans and other raw ingredients through the chocolate making process and including the final chocolate product.

The Unity Chocolate Analyzer contains calibrations for milk chocolate, dark chocolate, cocoa liquor, cocoa powder, cake, cocoa beans, nibs, and cocoa butter. Constituents predicted include fat, moisture, sugar, fatty acids and other constituents. Calibrations for ingredients such as milk powder are also available. All calibrations include 1 year of free calibration support.

Chocolate Analyzer

Robust and Ready to Use for At-Line or Lab Use

The Unity Chocolate Analyzer is robust enough to be used both at-line or in the lab. The system features a long life lamp and no air filter or fan, providing for very low cost of ownership. The built in computer is fully networkable and LIMS compatible, allowing for plant management to immediately access results. The Chocolate Analyzer is a valuable tool to control every step in the chocolate manufacturing process, from incoming ingredients to final product verification.

The Chocolate Analyzer package includes the following:

- SpectraStar 1400 XT (1400 nm - 2600 nm) with built in 17" touch screen, prediction software, multi-cup adapter, US-RTIR-0016 adapter ring, Unity TAS Standards
- 5 Ring cups US-ISIH-0007 for the analysis of powders
- Large cup US-LGOP-0001 for analysis of cocoa beans
- Small cup US-SRCP-0025 for analysis of powders and cocoa butter
- Gold Reflector US-TSTD-0005 for analysis of cocoa butter and liquor
- Unity Chocolate Calibration package for cocoa, chocolate and ingredients



HOW TO ORDER


The Chocolate Analyzer package includes everything you need to start analyzing cocoa and chocolate right out of the box.

CHOCOLATE ANALYZER

US-14XT-CHOC

SpectraStar 1400 XT

(1400 nm – 2600 nm)

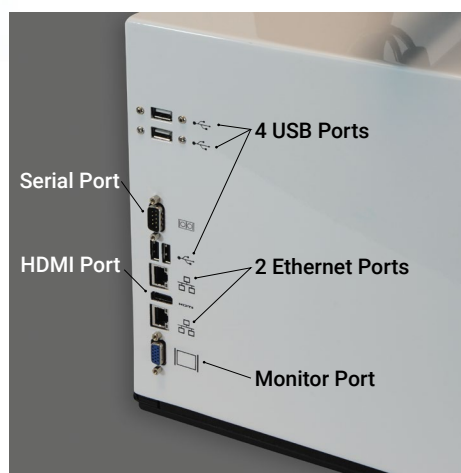

Specifications Table

Technology	Pre-dispersive scanning monochromator with nominal bandwidth of 10 nm (FWHM)
Wavelength Range	1400 - 2600 nm
Light Source	Tungsten halogen lamp with MTBF rating of 10,000 hours (User changeable)
Detector	Custom dual cooled, extended InGaAs detectors (2)
Analysis Time	10 - 30 seconds
Data Interval	1.0 nm
Photometric Noise	< 0.02 mAu 1400 - 2600 nm

User Interface

Built-in Computer	Windows® 7 Embedded, Solid State HDD, 17" high resolution touch screen monitor
Software	UScan™ prediction and data management package
Network Capabilities	OPC Compliant; LIMS Compatible; 5 USB ports, 2 Ethernet (RJ45)

Dimensions and Power Requirements

Size	330 x 381 x 508 mm (with display) 330 x 381 x 270 mm (without display)
Weight	18 kg, 40 lbs.; 22 kg, 48 lbs. with monitor
Voltage	100-240 V AC *, frequency 50-60 Hz, Class 1, protective earth

Unity Scientific Asia Pacific Pty Ltd
Unit 7/ 84-90 Old Bathurst Rd, Emu Plains. NSW, Australia
Ph: 02 4735 1813
E: sydney@unityscientific.com.au
www.unityscientific.com.au

Unity Scientific, InfoStar, SpectraStar XL UScan, and UCal are all trademarks of Westco Scientific Instruments, Inc., d/b/a Unity Scientific. All other trademarks are the property of their respective companies. 0901B ©Copyright 2016 All rights reserved. Form #ES_004_05.20.16